

Quant-iT™ RNA Assay Kit, Broad Range (Q10213)

Protocol summary

1. Equilibrate assay components to room temperature.
2. Make the working solution by diluting Quant-iT™ RNA BR reagent 1:200 in Quant-iT™ RNA BR buffer.
3. Load 200 µl of the working solution in each microplate well.
4. Add 10 µl of each of the Quant-iT™ RNA BR standards to separate wells, and mix well.
5. Add 1–20 µl of each unknown RNA sample to separate wells, and mix well.
6. Measure fluorescence using a microplate reader (excitation/emission maxima ~644/673 nm).
7. Use a standard curve to determine RNA amounts.



www.invitrogen.com